

PRECELLYS® 24 TOUCH



Intuitive Tissue Homogenizer

- **Easy to use:** touchscreen and exclusive tube locking system
- **Efficient:** powerful 3D motion, up to 6,800 rpm
- **Reproducible:** validated protocols available online
- **High quality:** robust, ergonomic & safe

PRECELLYS® 24 TOUCH

Intuitive Homogenizer




Precellys® 24 Touch is the ideal lab homogenizer to prepare any kind of biological sample, from the softest to the hardest, regardless of its origin (animal tissues, plant tissues, microorganisms...). It can treat up to 24 samples at a time and has been validated on more than 1,500 different samples: from brain, muscle and heart samples, to very hard material such as bones, hair and seeds.

The intuitive touchscreen interface of Precellys® 24 Touch makes it really easy to use. Thanks to its exclusive vacuum-based tube locking system and its specific design, Precellys® 24 Touch is compliant with safety requirements of the harshest lab environments, such as BSL3.

With its powerful figure-8 multi-directional motion, Precellys® 24 Touch provides high energy to the beads contained in the tubes, allowing the homogenization of the samples in a few seconds (~30 sec) at speeds ranging from 4000 to 6800 rpm. Samples can then be processed for further analysis amongst a wide variety of techniques dedicated to DNA/RNA (PCR, electrophoresis, NGS, chips, etc), protein (blotting, mass spectrometry, etc.), or metabolites analyses (chromatography, etc.).

Applications

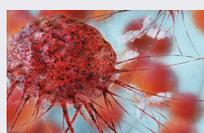

Cancer research



Agronomy



Forensic



Pharmacology




Environment




Microbiology


Precellys® 24 Touch - Quick User Mode

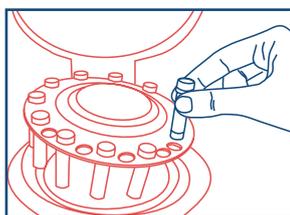


1. Load from one to twenty-four 0,5 or 2ml tubes

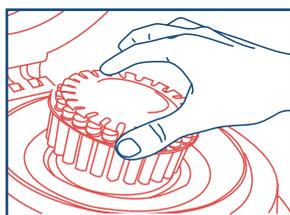

2. Position the patented system for quick and safe locking

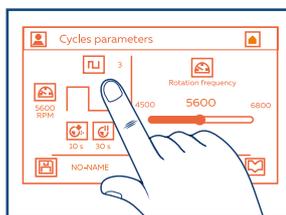


3. Set speed (4,000 - 6,800 rpm), time and cycle

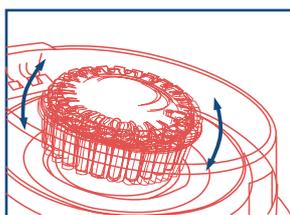

4. 3D bead-beating: homogenization is achieved in few seconds



5. The same efficiency in each tube enables the liberation of cell components

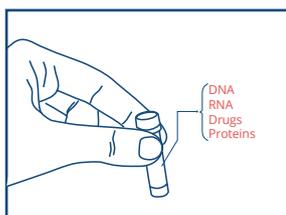

6. In 30 seconds, your sample is ready for analysis!



Technical specifications

CAPACITY:

1 to 24 tubes of 2 mL
1 to 24 tubes of 0.5 mL

SPEED RANGE:

4,000 rpm to 6,800 rpm
(by increments of 100 rpm)

PROGRAMS: Unlimited

NUMBER OF CYCLES: up to 10

TIME OF CYCLE: up to 90 sec

PAUSE BETWEEN CYCLES: 120 sec

SIZE:

L: 287 mm, W: 408 mm, H: 390 mm
(515mm with opened lid)

WEIGHT: 23 kg

NORMS: CE, FCC PART 15

ELECTRICITY:

90 - 250VAC / 50-60 Hz
Power: 1 kVA

NOISE: < 70dB

CATALOG NUMBER:

P002391-P24T0-A.0

Precellys homogenizers range

Discover the full range of products



Minilyt
for occasional use



Precellys® Evolution
the high-end versatile
homogenizer



Cryolys® Evolution
for sensitive molecules
(compatible with Precellys® Evolution)



**Lysing
kits consumables**